

## ProScope HR Quick Start

### Macintosh:

1. Make sure you have Mac OS 10.4.8 and Quicktime 7 or later installed on your computer.\*
2. Launch (double-click) the ProScope HR software installer found on the included CD-ROM. Macintosh folder>ProScope HR Installer.
3. Please quit iChat and unplug (or turn off) any external video camera before plugging in the ProScope HR USB microscope.
4. Plug the ProScope HR USB microscope into a factory installed USB 2.0 or USB 1.1 port on your computer.\*\*
5. Launch (double-click) the ProScope HR software application located in your applications folder. Now you can begin using your ProScope HR USB microscope.\*\*\*

### Windows:

1. Make sure you have Windows XP Service Pack 2 and QuickTime 7 or later installed on your computer.\*
2. Launch (double-click) the ProScope HR software installer located on the included CD-ROM. Windows folder>InstallProScopeHR.
3. Plug the ProScope HR USB microscope into a USB 2.0 or USB 1.1 port on your computer.\*\*
4. Launch (double-click) the ProScope HR software application from your desktop ProScope HR icon or "Programs" Bodelin folder. Now you can begin using the ProScope HR USB microscope.\*\*\*

\*QuickTime is included on the CD-ROM for your convenience. The Windows ProScope HR installer will automatically install QuickTime on your computer if necessary.

\*\*USB 2.0 port recommended. A USB 1.1 port may be used but only 640x480 (VGA) and 320x240 (QVGA) modes are supported and at a much reduced frame rate. Refer to the ProScope HR technical specifications.

\*\*\*Make sure you plug in the ProScope HR USB microscope BEFORE launching the ProScope HR software application.

Important READ ME FIRST files and additional documentation are available on the CD-ROM provided or online at [www.proscopehr.com](http://www.proscopehr.com).